

Instructions for the most professional way to use FlexiQuot labels

1Cryobio produces the labels for his customers with the greatest possible care. The quality of the product must also be safe at the costumers place during the processing. Therefore 1Cryobio recommends the handling of all label according to the following parameters:

Storage:


Keep the rolls in the original packaging until to processing and store then around 20° C ± 2°C and 50% relative humidity in a well-aired dark room.

Processing:

- Place rolls carefully on mandrel, avoid bumping into it.
- Ensure a continuous and jerk-free unwind tension
- Avoid tightening pulls within the rolls
- Use the rolls in one processing step without longer downtimes
- Use a detour rollers of greatest possible diameter
- Ensure a near traction precedence
- Limit the use of labels to one single workflow. In case of multiple uses the rewind of the roll can lead problems in the process flow.
- Use only course guidance at donor machines in the original state

Durability / Processing time frame:

The roll shall be processed best within 12 months from the production date which is indicated on the label in the reel. Within this period 1Cryobio is able to guarantee the requested manufacturing quality.

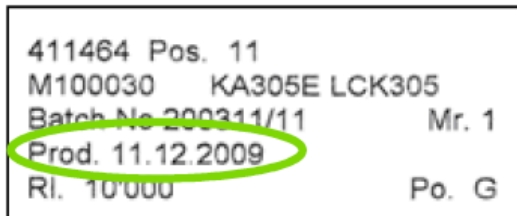

Warranty

1Cryobio rejects any warranties or liability for improper treatment of its label rolls.

1Cryobio AG